


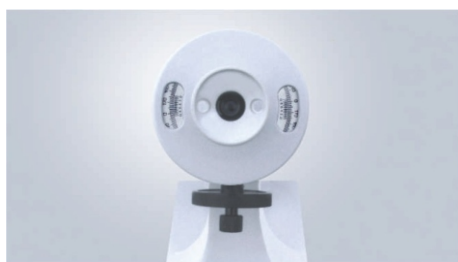

Rotate the vernier knob till the visual field to equal brightness. (Zero point)



**Full-circle manual polarimeter** is ideal for determining concentration levels of samples for application in the medical, educational, pharmaceutical, and cosmetic industries etc.

#### Easy Measurement

The polarimeter have 589.44nm sodium lamp. After powered on, the polarized light, beam is immediately radiate to the filter, polarizer etc. You can observe distinct visual fields through eyepiece of instrument. Put the sample tube into measuring field chamber. Rotate the vernier knob again, you can view equal brightness. Read the scale from magnifier that is measured value.

#### Specific optical Rotation Measurement

Specific optical rotation is constant of substance, you can measure it by 100mm glass tube when temperature of solution is at 20°C

#### Purity Measurement

If you want to know sample's purity you need only measure it's the angle of rotation and calculate it.

#### Concentration Measurement

For substance of known specific optical rotation, you can measure the angle of rotation and calculation concentration accordingly.

#### International Sugar Scale

According to international sugar scale, the 26 grams of pure sugar in solution, the angle of rotation of solution is 34.262°. By rule, you can calculate the purity of sugar.

| Model              | APM - 4                 |
|--------------------|-------------------------|
| Measurement Range  | -180° to +180°          |
| Scale Value        | 1°                      |
| Vernier            | ±0.05°                  |
| Magnifier          | 3X                      |
| Light Source       | Sodium Lamp             |
| Wavelength         | 589.44nm                |
| Parameter          | Manual                  |
| Round Cuvetter     | 100 or 200mm            |
| Stable Time        | 10 minutes              |
| Range of Accuracy  | 0 ~ ±180° (±0.05°)      |
| Power Requirements | 230 Vac, ± 20%, 50-60Hz |
| Dimension (LxWxH)  | 500 x 135 x 330mm       |
| Weight Approx.     | 6.5 kg                  |

WORLDWIDEMARKETING

**ACZET PVT. LTD.**

Unit E2, Plot No. 15, WICEL Estate, Opp Seepz Gate Nol. 1, Andheri (E), Mumbai 400 093. (INDIA)  
 Tel. : +91-22-4243 7700 • Fax : +91-22-4243 7800 • E-mail : sales@aczet.com • Web : www.aczet.com